


LectroPol-5


 **Struers**

Automatic, micro-processor controlled electrolytic polishing and etching of metallographic specimens



- Scanning function for easy determination of parameters
- Built-in safety features
- Database with methods for various materials
- Short polishing times and maximum reproducibility

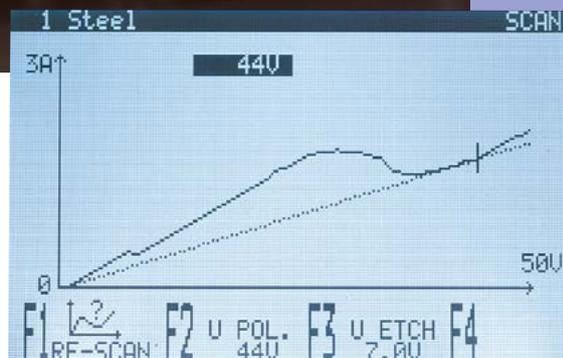

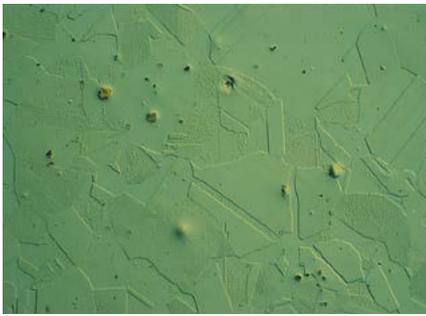

Stainless Steel,
500x, DIC




Copper,
500x, DIC

Scanning function for easy determination of parameters

LectroPol-5 is equipped with a unique scanning function. The sample is placed on the polishing table and a pre-defined voltage range is scanned to determine the current density curve.

This curve is used to define the correct voltage for both polishing and etching.

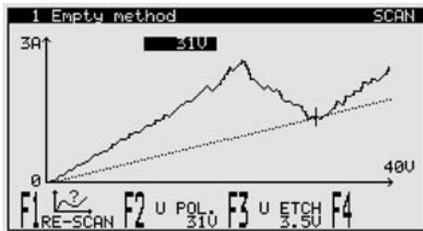

Instead of time consuming trial and error testing of parameters it is now easy to find the correct settings after a single scan of a new sample material.

Electrolytic polishing has now lost a lot of its "mystery" and is used in a more scientific way to prepare metallographic samples.

Built-in safety features

With LectroPol-5, all functions are controlled through the advanced software. An increase in electrolyte temperature above a pre-defined limit will automatically result in either a warning or shut-down of the procedure in progress.

Also voltage and current are monitored constantly and if the limits of Lectro-Pol-5 are reached, settings are reduced automatically.

Thus maximum safety for both operator and equipment is provided.

Database with methods for various materials

10 polishing/etching methods for various materials are included with LectroPol-5. A whole range of materials can thus be prepared immediately, without any lengthy and time consuming trials. These methods can also be used as starting point for the development of methods for other materials. 20 user-defined methods can also be saved in the database of LectroPol-5.


Short polishing times and maximum reproducibility

Electrolytic polishing and etching have the advantage of very short preparation times compared to normal mechanical specimen preparation. Microprocessor control and database functions result in always the exact same parameters being used. LectroPol-5 combines both short preparation times and high reproducibility and is thus the ideal choice for fast quality control requirements.

LectroPol-5 consists of two separate units, the control unit and the polishing unit.

The control unit

Both the power supply and the programming and monitoring functions are incorporated in the control unit. The unit is operated on the touch pad and on the backlit graphic display all parameters are presented. Upon pressing the Start button, the display

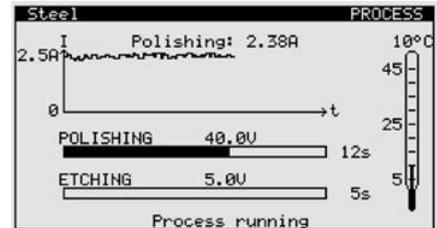


changes, and the ongoing process can be monitored. Parameters like current, electrolyte temperature and elapsed polishing/etching time are displayed. Both the display of parameters before and the display of actual values after process start provide an exceptional amount of immediately useable information. Deviations from one process to the next can be detected straightaway, and necessary measures can be taken. Required changing of electrolyte can be detected before changes of the polishing result will be noticed.


The polishing unit

The design as a separate unit makes it possible to place the polishing unit away from the control unit. If required the polishing unit can be set up in a fume cabinet while the control unit is set up outside. With an additional start and stop key the process can also be started directly from the polishing unit. The electrolyte is stored in easily exchangeable electrolyte containers which are inserted into the polishing unit.

The pump, driven like a magnetic stirrer, and the cooling coils are lowered into the container. On the polishing table the sample is positioned and electric contact is established through the anode arm. Two different polishing units are available.

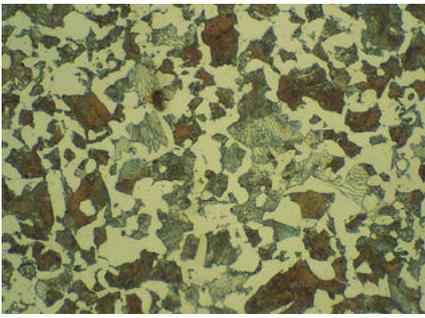

Medium Carbon Steel, 500x

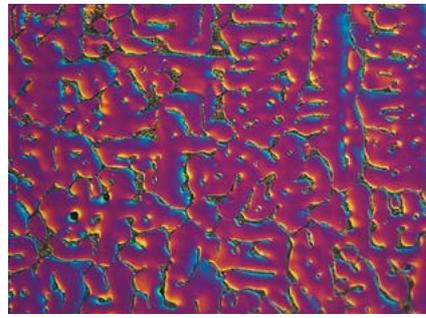

Aluminium alloy cast, 200x, DIC

With advanced cooling control

LectroPol-5 is equipped with a built-in measuring and control system, constantly monitoring the electrolyte temperature. As soon as a pre-set temperature is reached, a solenoid valve in the polishing table is activated, allowing water from the water mains or liquid from an external cooling unit to run through the cooling coil in the electrolyte container. This saves precious resources as the flow of cooling water is limited to the periods where the electrolyte temperature is above a pre-set limit.

If no water or cooling unit is connected, or the electrolyte temperature for some other reasons increases further, a warning can be displayed and eventually LectroPol-5 will shut off the process automatically.

For polishing at sub-zero temperatures

With this unit, the stainless steel cooling coils are led directly to the back of the polishing unit. There they can be connected to an external cooling unit. The external cooling unit takes over the temperature control and regulation of the electrolyte. With this set-up, materials can be polished at sub-zero temperatures.

External etching

LectroPol-5 is also prepared for external etching. This is often used when different electrolytes have to be used

for polishing and etching. Also with external etching a time can be set, and LectroPol-5 will automatically detect when the specimen is lowered into the external etching bowl. Then the countdown is running, and the voltage is terminated when the pre-set time has expired.

Accessories

Apart from the standard equipment, we can offer additional accessories for special applications.

Switchbox


To avoid the constant cleaning when changing between different electrolytes, LectroPol-5 can be equipped with two polishing tables and a switchbox to change between these two tables.

Extension arm

For the preparation of high specimens an extension arm can be mounted.



The transparent masks facilitate monitoring the electrolyte level and the adjustment of the electrolyte flow.

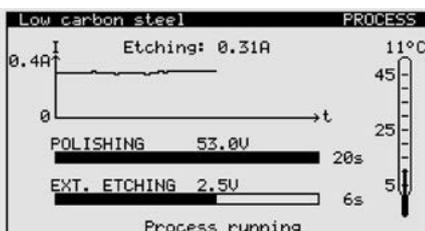

Technical Data

Power Supply

50/60Hz (max. load 9.8A)	1 x 100V / 120V
50/60Hz (max. load 4.9A)	1 x 220V / 240V

Output voltage/current

Polishing	0 - 100V	(1V steps) / 6A
Etching	0 - 25V	(0.5 V steps) / 6A
External Etching	0 - 15V	(0.5 V steps) / 1.5A

Software and Electronics

Display	128 x 240 dots (16 x 40 characters)	
Controls	Touch pad	
Database	10 Struers methods + 20 user definable methods	

Dimensions and Weight

	Control unit	Polishing unit
Width	385 mm / 15.2"	220 mm / 8.7"
Depth	350 mm / 13.8"	350 mm / 13.8"
Height	160 mm / 6.3"	160 mm / 6.3"
Weight	18 kg / 40 lbs	6.3 kg / 14 lbs

Specifications

LectroPol-5

Fully automatic, microprocessor controlled electrolytic polishing and etching apparatus. Complete with control unit, polishing unit (04936201) and external etching unit (02250121). One set of masks (04936915) and two electrolyte containers (04936902) included.

Control unit with external etching unit

Polishing unit with one set of masks (04936915) and two electrolyte containers (04936902)

Polishing unit for low temperature, with one set of masks (04936915), one cathode (04936908) and two electrolyte containers (04936902). External cooling unit and tubing for connection to the external cooling unit are not included and have to be sourced alternatively. Internal dia. of tubing to be used, 6 mm

Switchbox for the connection of two polishing units to a LectroPol-5 Control Unit.

Extension Arm, for the preparation of specimens up to 100 mm high. Can be mounted on pol. unit 04936202 and on pol. unit 04936201 from serial no. 4935250

1 set of masks (0,5, 1, 2, 5 cm² and 1 without aperture.)

1 set of 5 masks 0,5 cm²

1 set of 5 masks 1 cm²

1 set of 5 masks 2 cm²

1 set of 5 masks 5 cm²

1 set of 5 masks without aperture

Electrolyte container with lid

External etching unit

Cat. no.

04936133

04936233

04936201

04936202

04936911

04936909

04936915

04936916

04936917

04936918

04936919

04936920

04936902

02250121

List of Electrolytes

A2	A3	D2
Aluminium	Martensitic Steel	Brass
Iron	Molybdenum	Copper
Nickel	Stainless Steel	Gold
Silver	Titanium	
Stainless Steel	Vanadium	
Steels	Zirconium	
Tin		
Titanium		

Struers' equipment is in conformity with the provisions of the applicable International Directives and their appurtenant Standards. (Please contact your local supplier for details)

Struers' products are subject to constant product development. Therefore, we reserve ourselves the right to introduce changes in our products without notice.

AUSTRALIA & NEW ZEALAND

Struers Australia
 27 Mayneview Street
 Milton QLD 4064
 Australia
 Phone +61 7 3512 9600
 Fax +61 7 3369 8200
 info.au@struers.dk

BELGIQUE (Wallonie)

Struers S.A.S.
 370, rue du Marché Rollay
 F- 94507 Champigny
 sur Marne Cedex
 Téléphone +33 1 5509 1430
 Télécopie +33 1 5509 1449
 struers@struers.fr

BELGIUM (Flanders)

Struers GmbH Nederland
 Zomerdijk 34 A
 3143 CT Maassluis
 Telefoon +31 (10) 599 7209
 Fax +31 (10) 597201
 netherlands@struers.de

CANADA

Struers Ltd.
 7275 West Credit Avenue
 Mississauga, Ontario L5N 5M9
 Phone +1 905-814-8855
 Fax +1 905-814-1440
 info@struers.com

CHINA

Struers Ltd.
 No. 1696 Zhang Heng Road
 Zhang Jiang Hi-Tech Park
 Shanghai 201203, P.R. China
 Phone +86 (21) 6035 3900
 Fax +86 (21) 6035 3999
 struers@struers.cn

CZECH REPUBLIC & SLOVAKIA

Struers GmbH Organizační složka
 vědeckotechnický park
 Přílepská 1920,
 CZ-252 63 Roztoky u Prahy
 Phone +420 233 312 625
 Fax +420 233 312 640
 czechrepublic@struers.de
 slovakia@struers.de

DEUTSCHLAND

Struers GmbH
 Carl-Friedrich-Benz-Straße 5
 D-47877 Willich
 Telefon +49 (0) 2154 486-0
 Fax +49 (0) 2154 486-222
 verkauf@struers.de

FRANCE

Struers S.A.S.
 370, rue du Marché Rollay
 F-94507 Champigny
 sur Marne Cedex
 Téléphone +33 1 5509 1430
 Télécopie +33 1 5509 1449
 struers@struers.fr

HUNGARY

Struers GmbH
 Magyarországi Fióktelep
 Tatai ut 53
 2821 Gyermely
 Phone +36 (34) 880546
 Fax +36 (34) 880547
 hungary@struers.de

IRELAND

Struers Ltd.
 Unit 11 Evolution@ AMP
 Whittle Way, Catcliffe
 Rotherham S60 5BL
 Tel. +44 0845 604 6664
 Fax +44 0845 604 6651
 info@struers.co.uk

ITALY

Struers Italia
 Via Monte Grappa 80/4
 20020 Arese (MI)
 Tel. +39-02/38236281
 Fax +39-02/38236274
 struers.it@struers.it

JAPAN

Muramoto Struers K.K.
 Takara 3rd Building
 18-6, Higashi Ueno 1-chome
 Taito-ku, Tokyo 110-0015
 Phone +81 3 5688 2914
 Fax +81 3 5688 2927
 struers@struers.co.jp

NETHERLANDS

Struers GmbH Nederland
 Zomerdijk 34 A
 3143 CT Maassluis
 Telefoon +31 (10) 599 7209
 Fax +31 (10) 5997201
 netherlands@struers.de

NORWAY

Struers ApS, Norge
 Sjøskogenveien 44C
 1407 Vinterbro
 Telefon +47 970 94 285
 info@struers.no

ÖSTERREICH

Struers GmbH
 Zweigniederlassung Österreich
 Betriebsgebiet Puch Nord 8
 5412 Puch
 Telefon +43 6245 70567
 Fax +43 6245 70567-78
 austria@struers.de

POLAND

Struers Sp. z o.o.
 Oddział w Polsce
 ul. Jasnogórska 44
 31-358 Kraków
 Phone +48 12 661 20 60
 Fax +48 12 626 01 46
 poland@struers.de

ROMANIA

Struers GmbH, Sucursala Bucuresti
 Str. Preciziei nr. 6R
 062203 sector 6, Bucuresti
 Phone +40 (31) 101 9548
 Fax +40 (31) 101 9549
 romania@struers.de

SCHWEIZ

Struers GmbH
 Zweigniederlassung Schweiz
 Weissenbrunnstraße 41
 CH-8903 Birmensdorf
 Telefon +41 44 777 63 07
 Fax +41 44 777 63 09
 switzerland@struers.de

SINGAPORE

Struers Singapore
 627A Aljunied Road,
 #07-08 BizTech Centre
 Singapore 389842
 Phone +65 6299 2268
 Fax +65 6299 2661
 struers.sg@struers.dk

SPAIN

Struers España
 Camino Cerro de los Gamos 1
 Building 1 - Pozuelo de Alarcón
 CP 28224 Madrid
 Teléfono +34 917 901 204
 Fax +34 917 901 112
 struers.es@struers.es

SUOMI

Struers ApS, Suomi
 Hietalahdenranta 13
 00180 Helsinki
 Puhelin +358 (0)207 919 430
 Faksi +358 (0)207 919 431
 finland@struers.fi

SWEDEN

Struers Sverige
 Box 20038
 161 02 Bromma
 Telefon +46 (0)8 447 53 90
 Telefax +46 (0)8 447 53 99
 info@struers.se

UNITED KINGDOM

Struers Ltd.
 Unit 11 Evolution @ AMP
 Whittle Way, Catcliffe
 Rotherham S60 5BL
 Tel. +44 0845 604 6664
 Fax +44 0845 604 6651
 info@struers.co.uk

USA

Struers Inc.
 24766 Detroit Road
 Westlake, OH 44145-1598
 Phone +1 440 871 0071
 Fax +1 440 871 8188
 info@struers.com